

X-Max

The largest area SDD

SIZE

IT MATTERS

ALL WALLS?

20mm²

50mm²

80mm²



For **accuracy** / For **nano-analysis** / For **throughput** / For **productivity**

The Business of Science[®]




X-Max

Large area SDD

Analysis with a large area analytical SDD benefits from:

- Simultaneous imaging and analysis without compromise
- Large active area with superb analytical resolution
- Practical nanoanalysis at productive count rates
 - High count rates, even at low kV
 - High count rates, even at small spot sizes
- The correct results

A traditional 10mm² detector has a small capture angle = fewer counts
A modern large area detector has a large capture angle = many counts

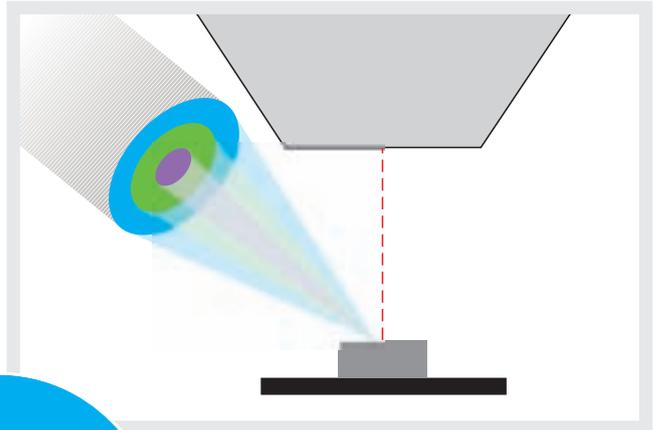


20mm²

50mm²

80mm²

X-Max delivers:

- Up to 1,000% greater productivity
- Real 200,000 cps analysis
- Usable nanoscale analysis
- MnK_α resolution down to 123eV - even at 80mm²

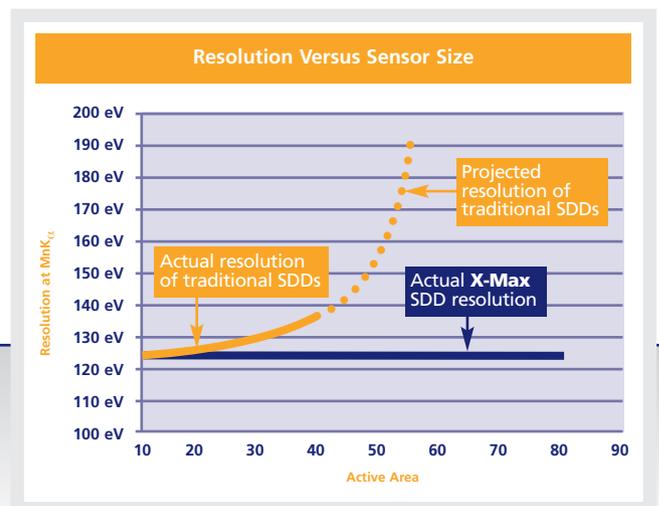

— Traditional SDD
— X-Max SDD

X-Max The largest area SDD

X-Max

Productivity and accuracy combined


20mm², 50mm², and 80mm² SDDs with identical world-beating analytical performance

Now you can have count rate, imaging, and analytical performance all at the same time:

- Up to 80mm² active area
- Count rates > 500,000 cps
- Throughput > 200,000 cps
- MnK_α typically 125eV

Maximum counts from larger detectors with increased solid angle:

- Count for less time - increase productivity
- Or count for the same time and get more precise results
- Or gather data at much lower beam currents meaning:
 - No sample damage
 - Longer filament tip lifetimes
 - Reduced contamination
 - Better spatial resolution
- Practical nanoanalysis with productive count rates at low accelerating voltages

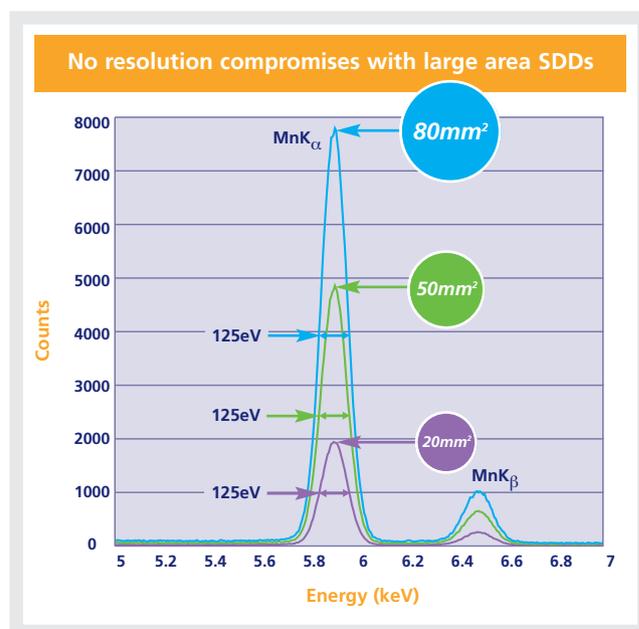

X-Max analytical SDDs offer the best solution for productive analysis

X-Max features


- Unique large area SDD sensors
- Discrete PentaFET[®] design - accuracy at all sizes
- Sensor enclosed in self-contained vacuum assembly - no oxygen X-ray absorption
- Unique electron trap for maximised solid angle
- Only one pulse processing channel required
- Tube diameter no bigger than a 10mm²

SIZE
IT MATTERS
ALL MATTERS?

X-Max

Large area SDD

Why size matters

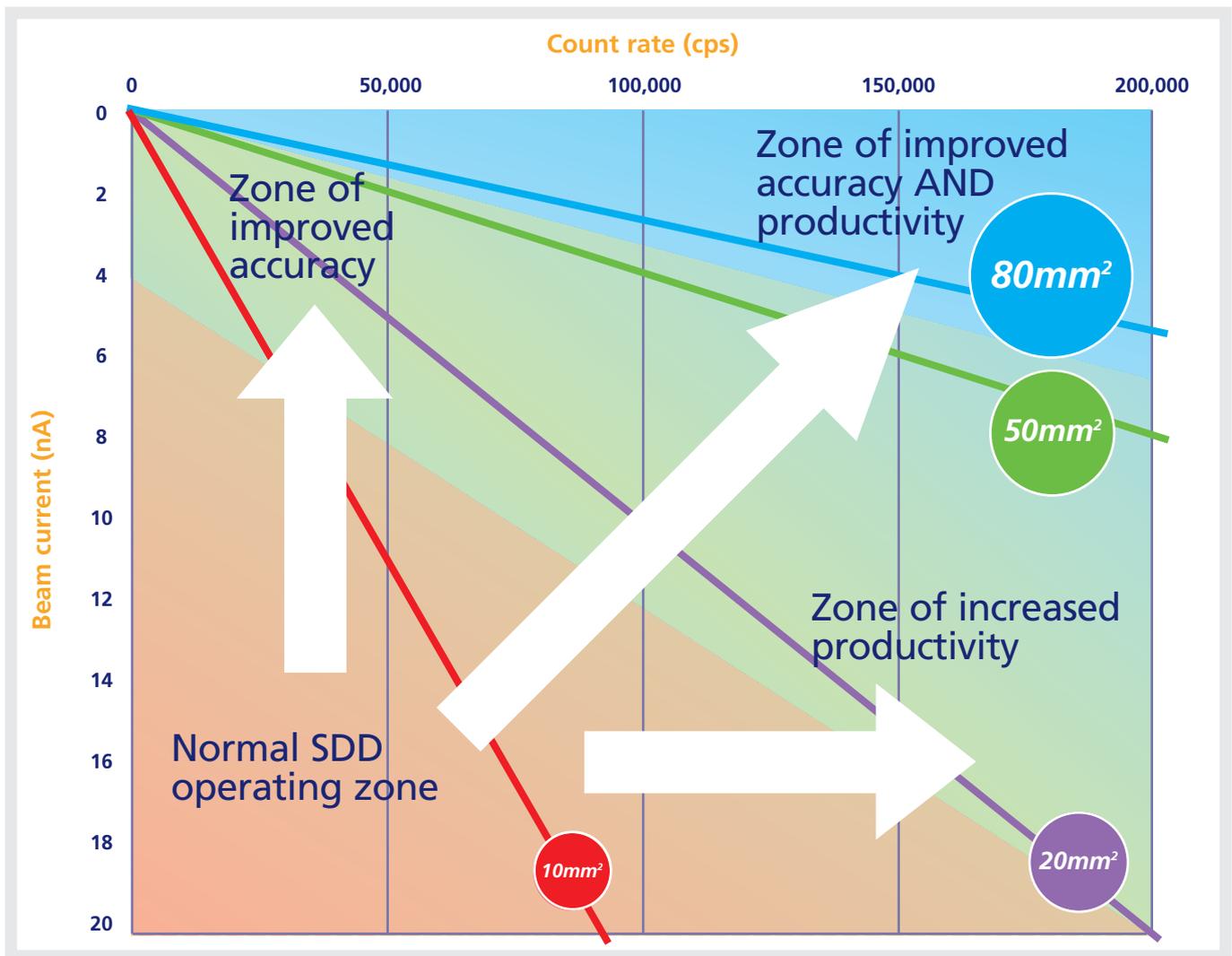

Discover the benefits of working with large SDDs.

Previously, users had to choose between:

1. Working at low beam currents to maintain accuracy
2. Working with high beam current to get high count rates, risking beam damage and contamination

Now X-Max offers users two more options:


3. Work at low beam currents with high count rates and improved accuracy
4. Or decrease analysis times

X-Max The largest area SDD

X-Max

Size and productivity


Benefit immediately with increased counts from **X-Max**

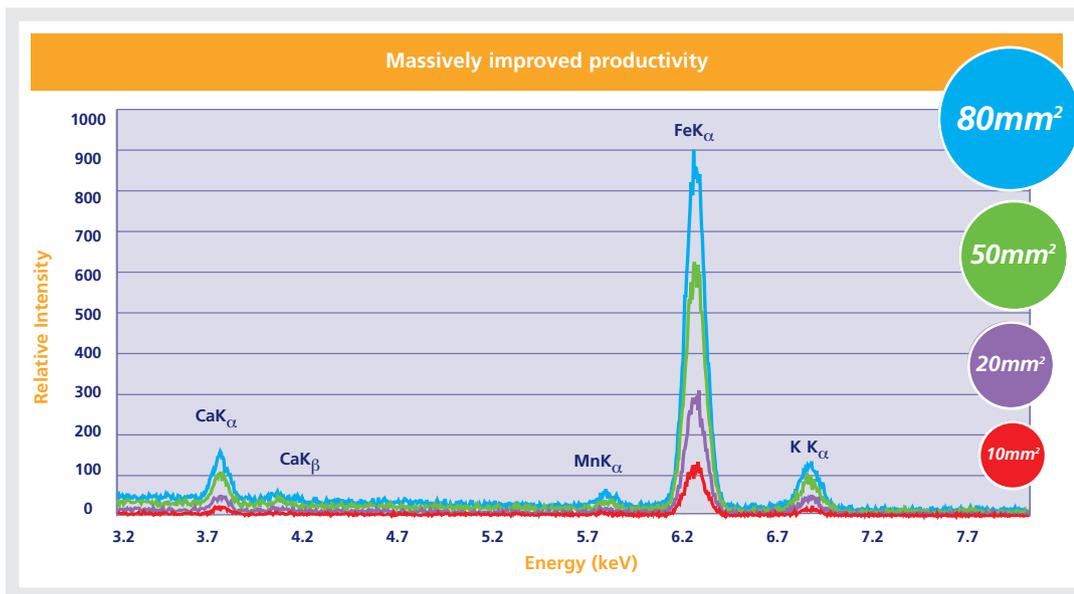

These almandine garnet spectra were collected using a 10mm² SDD detector and the new 20mm², 50mm², and 80mm² X-Max detectors. The larger X-Max detectors collect more X-rays under the same collection conditions.

True 100,000 cps performance

Many more analyses can be performed with 80mm² detectors at 100,000 cps

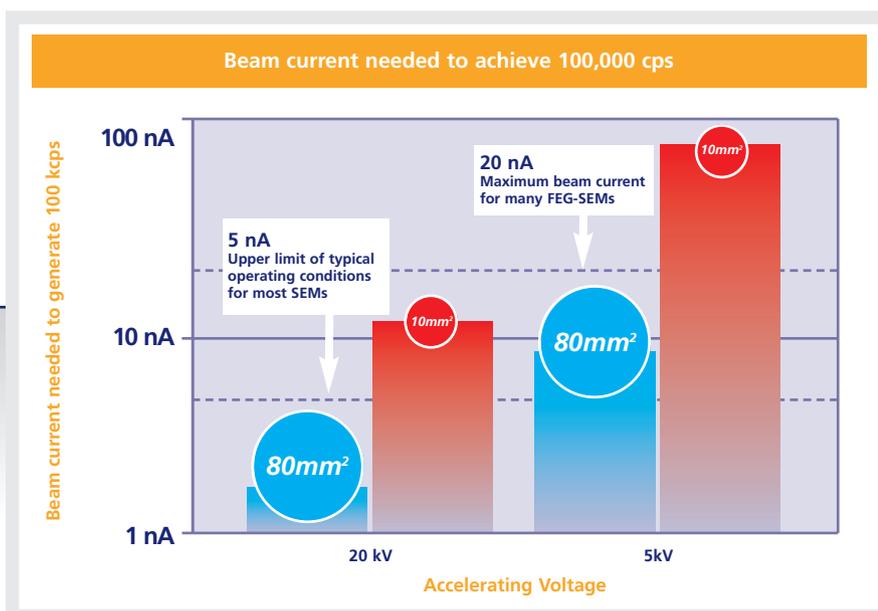

This logarithmic scale bar chart shows how high the beam current needs to be to achieve 100kcps on a traditional 10mm² SDD at 5kV and 20kV. It is a ten-fold increase over what is required to generate the same count rate on a new X-Max 80mm² SDD detector.

SIZE
IT MATTERS
ALL MATTERS?


X-Max

Large area SDD

The effect of beam current on analytical data

Data collected with standard 10mm² SDDs


Insufficient data for usable nanoanalysis

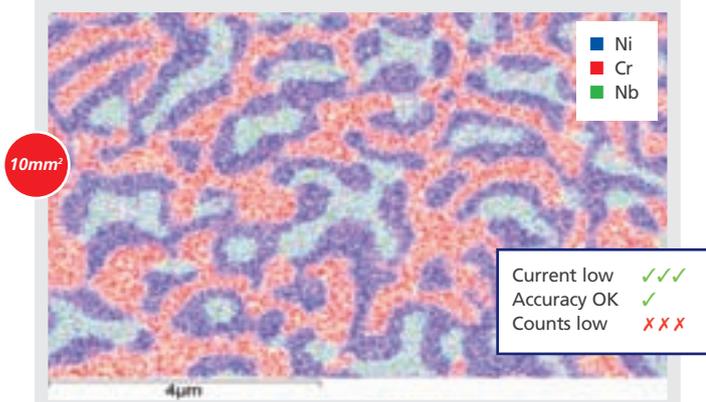

Some detail but very few counts

Data collected with a 10mm² detector for 45 minutes at 3.3nA. Insufficient data has been collected which means a poor map that is not usable for analysis.

High beam current reduces accuracy

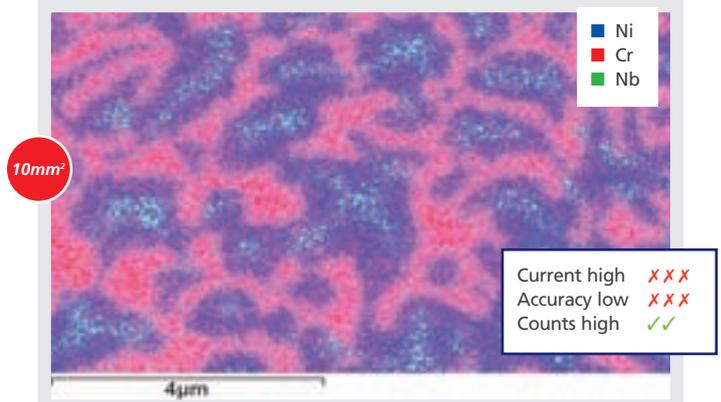


Lots of counts but no detail

Data collected with a 10mm² detector for 45 minutes at 36nA. The map shows lots of counts but the detailed information is blurred by a large spot size at this current.

Lots of counts and high quality data at low beam current

Data collected with 80mm² X-Max SDD

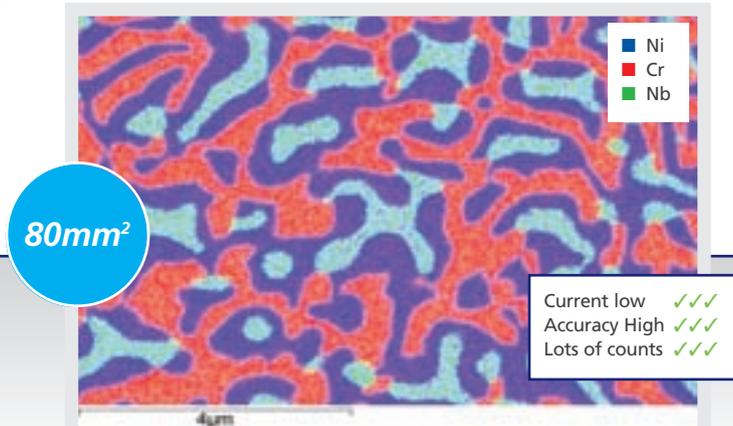

Lots of detail AND lots of counts


Data collected with an 80mm² detector for 45 minutes at 3.3nA. The mix map shows good spatial resolution and contains lots of data for good quantitative analysis.

X-Max

Size and accuracy

Increasing accuracy by lowering accelerating voltage

X-Max offers unparalleled benefits when working at low kV (and low count rates).

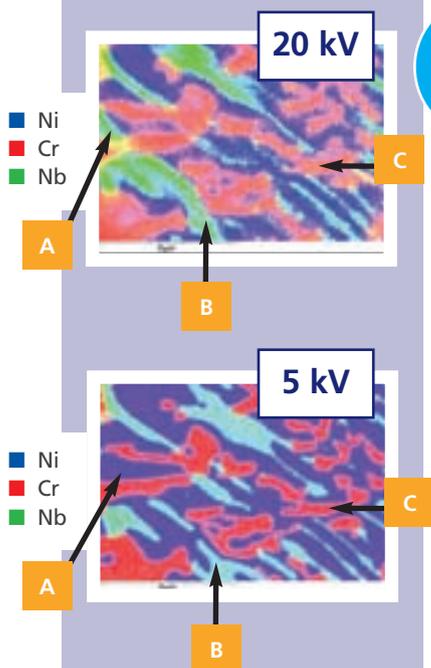

Maximise counts for accurate nanoanalysis with X-Max


These two maps were collected from the same area under the same conditions, changing only accelerating voltage. While superficially similar, points A, B, and C are actually very different. This is because reducing accelerating voltage (kV) reduces the interaction volume from which the X-rays are generated. As a result, the 20kV map shows X-ray signals from structures below the surface not seen in the 5kV map. Analysing at low kV benefits nanostructure analysis as it means that X-rays are only collected from the surface structures.

Larger detectors cause less surface damage

Beam damage on sensitive samples




This image is of a polymer sample, taken at 0.2nA. Features on the surface and striation from sample preparation can clearly be seen.

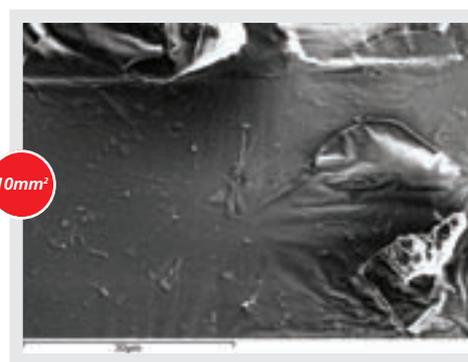

The same sample imaged at 2nA shows extensive surface damage and distortion as a result of the higher beam current.

Beam contamination

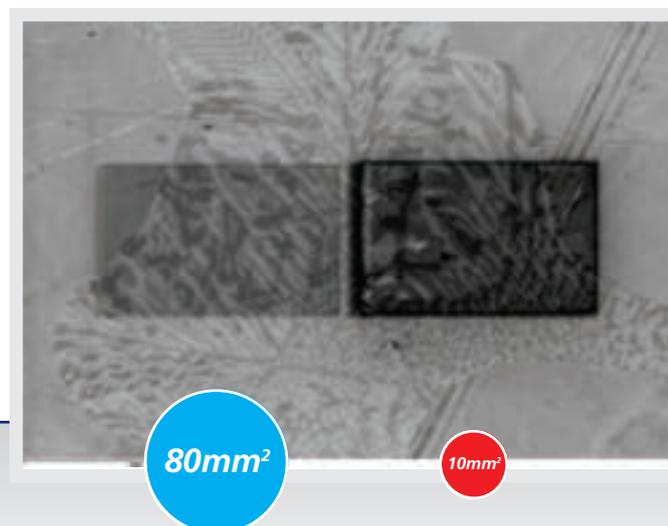

This image is of a sample from a nickel spinner bowl. Two areas have been examined, the one on the left using a large area detector, the one on the right with a traditional 10mm² detector. To collect the same data the sample had to be exposed for ten times longer with the smaller detector and as a result has far worse contamination.

X-Max

Technology leadership




Unique large area sensors

- Discrete PentaFET® design offers the following benefits:
 - Protects FET from X-ray damage*
 - Superior thermal design and FET gain stabilisation means detectors cool quickly and stabilise fast
 - Radial design for minimal ballistic deficit*
- Polycrystalline Si construction process for extreme low leakage



Unique magnetic trap design

- Allows large area sensor very close to samples
- Minimum sample to sensor distance = maximum capture angle
 - Typically 10 times greater than standard 10mm² SDDs
- Uses optimised rare earth element design

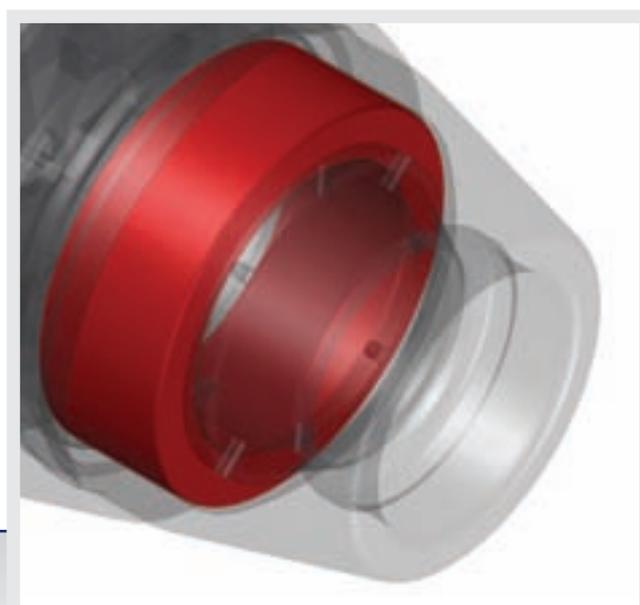

* For more information about silicon drift sensor technology and what makes an SDD detector analytical, go to www.oxinst.com/sdd to request a free copy of 'SDD Explained'

X-Max The largest area SDD


X-Max

Pushing the boundaries



Vacuum enclosed sensor

- No gas fill to absorb light elements
 - Nitrogen gas in other SDDs absorbs oxygen and other low energy X-rays
 - Vacuum enclosure increases analytical accuracy
- Long lifetime design
 - Uses OI vacuum technology
- Vacuum allows for greater cooling
 - Dramatically reduces thermal noise in large area SDDs
 - Not possible with gas-filled systems

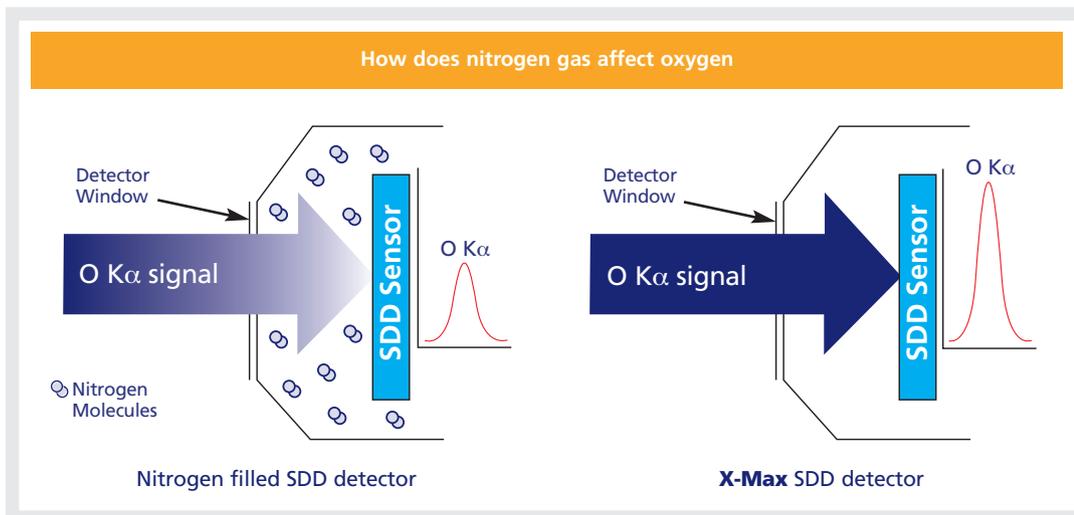

The oxygen X-rays are absorbed by the nitrogen molecules present in the 'vacuum' so only a fraction of the X-ray signal gets through to the sensor.

The oxygen X-rays are not absorbed as there is no nitrogen present so you can see the true shape and height of the oxygen peak.


On site 8-hour repair capability

- Includes window breakages

X-Max

Applications

Low energy mapping of nano-particles

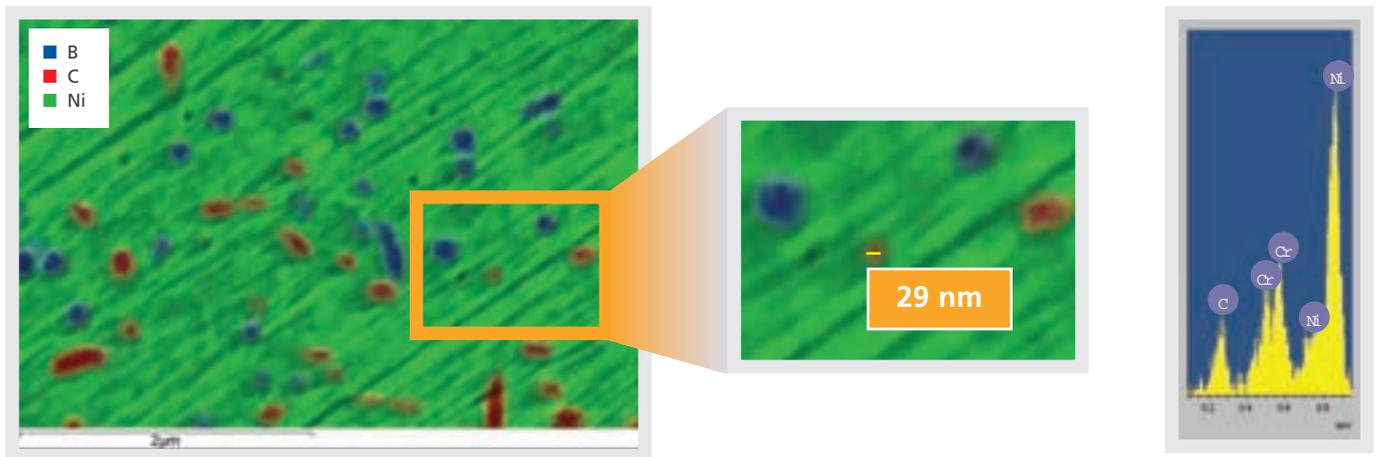


X-Max achieves excellent resolution. It also allows the user to do extraordinary nanoanalysis under conditions where other detectors struggle to produce sensible results. This means **X-Max** not only allows detection of small precipitates, but can map them as well.

Here we see a sample containing very small boride and carbide precipitates. The map clearly distinguishes between the different phases and even identifies the chemical composition of precipitates as small as 30nm. The spectrum above is from the highlighted 30nm carbide precipitate and includes data equivalent to a 0.5 second acquisition.

Very high count rate maps

Texture analysis of fracture surface of a granite sample. This map, taken at a count rate of 400,000cps with an **X-Max** 80mm² detector, shows clearly how it is possible to map the different chemical compositions of a rough surface to see how the textured surface and chemical composition of the surface relate to each other.

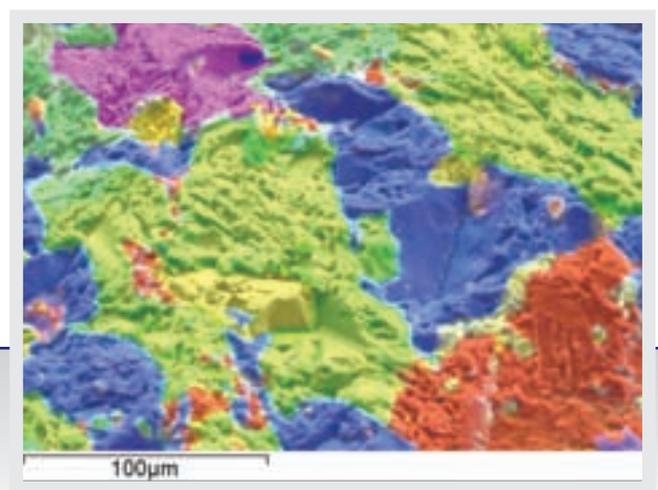


X-Max

Superb performance

Why does an **X-Max** EDS system give such exemplary performance?

- A great detector
 - Maximum real active area
 - Excellent resolution at high and low energies
 - Good throughput
 - ISO15632:2002 resolution compliance
- Excellent electronics
 - Highly linear analogue to digital conversion
 - Good rejection of simultaneous events (Pulse-pile-up)
 - Excellent signal to noise ratio
 - < 1eV change in peak resolution and position between 1 and 100,000cps
- Unique OI-owned interpretation algorithms
 - Complete - and correct - database of X-ray lines and edges
 - Software correction of sum-peaks
 - Highly accurate deconvolution, AutoID, and Quantitation

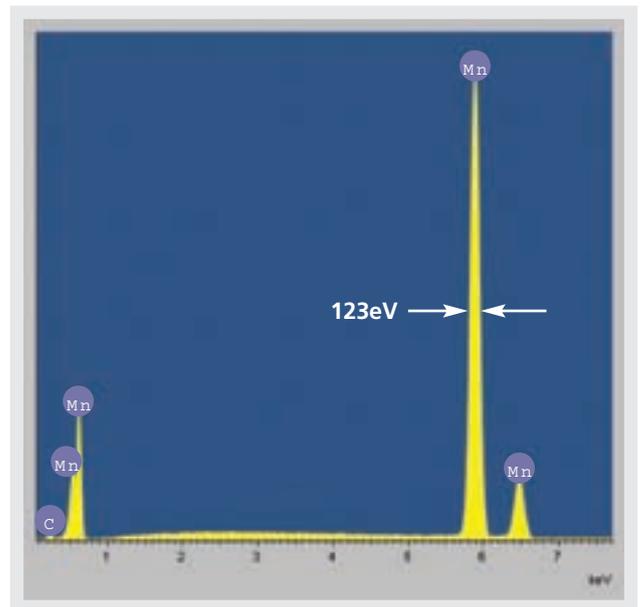

What the system doesn't have:

- Shelf, tail, and shift correction - only real data is analysed
- Fan cooling - no vibration
- Forbidden elements - all elements from Be to Pu can be analysed
- Identification of absent elements - no false positives

At 80mm², **X-Max** gives Mn_{K_α} resolution that other SDDs can't match. Here we see 123eV; 6 eV better than the guaranteed specification of 129eV.

(Spectrum courtesy of P Russell, Appalachian State University. Recorded at 20kV, 168pA).

X-Max

For **accuracy** / For **nano-analysis** / For **throughput** / For **productivity**

SIZE
IT MATTERS
ALL WALLS?

X-Max

In practice



Detector area	20mm ²	50mm ²	80mm ²
Resolution (eV):			
MnK _α resolution	129	129	129
Operating angle	0° to 45°		
Temperature range	10°C to 30°C		
Altitude	Sea level to 1,500m		

Comparison of detector type

	Si(Li)	Other SDD	X-Max
LN ₂ -free	NO	YES	YES
Maximum single sensor active area	30mm ²	30mm ²	80mm ²
Typical analysis conditions for most productive count rate			
Spot size	Small	Large	Small
Beam current	Low	High	Low
Accuracy at low kV	High	Low	High
Maximum throughput	Low	High	High
Productivity at low beam currents	Low	Low	High

Oxford Instruments, at High Wycombe, UK, operates Quality Management Systems approved to the requirements of BS EN ISO 9001. This publication is the copyright of Oxford Instruments Analytical Limited and provides outline information only which (unless agreed by the company in writing) may not be used, applied or reproduced for any purpose or form part of any order or contract or be regarded as a representation relating to the products or services concerned. Oxford Instruments' policy is one of continued improvement. The company reserves the right to alter, without notice, the specification, design or conditions of supply of any product or service. Oxford Instruments acknowledges all trade marks and registrations.

© Oxford Instruments Analytical Ltd, 2008. All rights reserved.



Part no: OIA/160/A/0908

As part of Oxford Instruments' environmental policy this brochure has been printed on FSC paper



Oxford Instruments NanoAnalysis

UK

Halifax Road, High Wycombe
Bucks, HP12 3SE England
Tel: +44 (0) 1494 442255
Fax: +44 (0) 1494 524129
Email: nanoanalysis@oxinst.com

China

Shanghai
Tel: +86 21 6360 8530/1/2/3
Fax: +86 21 6360 8535
Email: na.sales@oichina.cn

France

Saclay, Cedex
Tel: +33 (0) 1 69 85 25 24
Fax: +33 (0) 1 69 41 86 80
Email: analytical-info@oxford-instruments.fr

Germany

Wiesbaden
Tel: +49 6122 937 176
Fax: +49 6122 937 175
Email: nanoanalysis@oxford.de

Japan

Tokyo
Tel: +81 (0) 3 5245 3591
Fax: +81 (0) 3 5245 4466/4477
Email: nanoanalysis@oxinst.co.jp

Latin America

Concord, MA, USA
Tel: +1 978 369 9933
Fax +1 978 369 8287
Email: nanoanalysis@ma.oxinst.com

Singapore

Tel: +65 6337 6848
Fax: +65 6337 6286
Email: nanoanalysis@oxford-instruments.com.sg

Sweden

Lidingö
Tel: +46 8 544 81550
Fax: +46 8 544 81558
Email: nordiska@oxinst.com

USA - Oxford Instruments

Concord, MA, USA
Tel: +1 978 369 9933
Fax +1 978 369 8287
Email: nanoanalysis@ma.oxinst.com

www.oxford-instruments.com

